


# INFINITY STIRRER - MANUAL OPERATION

Short presses of the + and - buttons step through the following sequence.


From OFF two short - presses within two seconds will start the stirrer at 350 RPM.

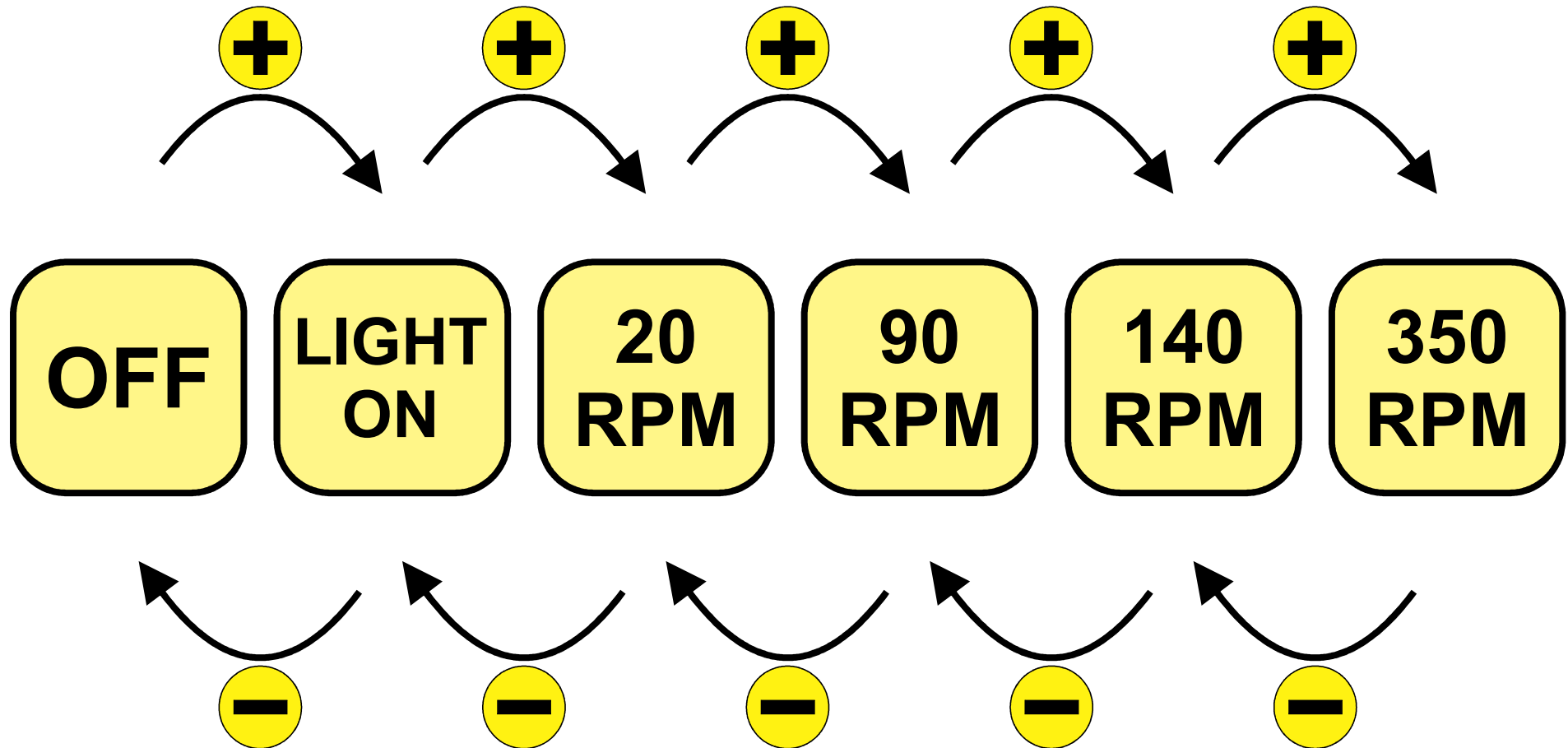

Pressing and HOLDING the + or - buttons will allow you to ramp the speed up or down to any value in the range of 20 RPM and 500 RPM.

For more information visit [www.boltac.co.nz](http://www.boltac.co.nz)

# INFINITY STIRRER - TABLET OPERATION

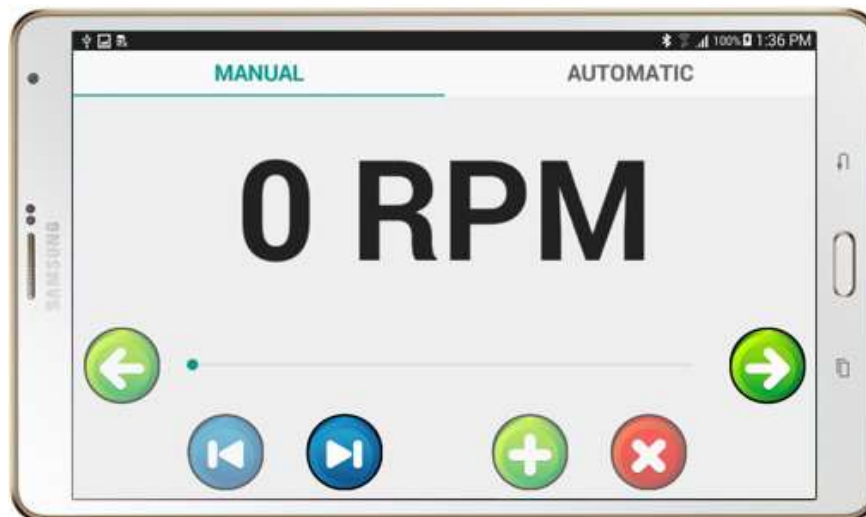

Your Infinity stirrer can be bluetooth controlled using an **ANDROID** tablet.  
To do this you will need to install the **BOLTAC STIRRER CONTROL** app from the google **PLAY STORE**.

Once installed you will need to **PAIR** your tablet with your stirrer using the **BLUETOOTH SETTINGS** in the **SETTINGS MENU**.

Look for a device named **INFINITY**. The pairing **PIN** is **1234**.

It is recommenced that you connect the **TABLET** to the **USB CHARGER** socket on the back of the stirrer control box while the stirrer is running.